

Compatible models for ORP60-DE replacement probe:

- **ORP60-Z**

How to replace the probe?

Screw off the probe ring, unplug the probe, plug in the new replacement probe (pay attention to the probe's position), and screw on the probe ring.

**Video tutorials are available at www.aperainst.de or on YouTube (Apera Instruments Europe Official Channel)

We warrant this electrode to be free from defects in material and workmanship and agree to repair or replace free of charge, at option of APERA INSTRUMENTS (Europe) GmbH any malfunctioned or damaged product attributable to responsibility of APERA INSTRUMENTS (Europe) GmbH for a period of six-month. This warranty does not apply to defects resulting from actions such as misuse (violation of the instructions in this manual or operations in the manner not specified in this manual), improper maintenance, or unauthorized repairs. Warranty period is the time limit to provide free service for the products purchased by customers, not the service life of the probe.

APERA INSTRUMENTS (Europe) GmbH
Wilhelm-Muthmann-Straße 18, 42329 Wuppertal, Germany
Tel. +49 202 51988998 | info@aperainst.de | www.aperainst.de



Il sensore di ricambio (elettrodo) APERA INSTRUMENTS ORP60-DE per la Serie 60-Z è compatibile con:

- **ORP60-Z**

- Rimuovere il cappuccio del sensore.
- Svitare l'anello sul sensore e allentare il sensore.
- Collegare il nuovo sensore al dispositivo. Assicurati che l'allineamento sia corretto.
- Riavvitare l'anello

** Puoi trovare video tutorial su www.aperainst.de o su YouTube (Apera Instruments Europe Official Channel).

APERA INSTRUMENTS (Europe) GmbH concede una garanzia di sei mesi su questo prodotto. La garanzia del produttore decorre dalla data del primo acquisto da parte del primo cliente finale (data fattura). La garanzia copre il perfetto funzionamento del prodotto. Se durante il periodo di garanzia dovessero emergere difetti del prodotto dovuti a errori di fabbricazione o di elaborazione, APERA INSTRUMENTS (Europe) GmbH riparerà il prodotto o la parte difettosa gratuitamente o (a sua discrezione) lo sostituirà. In particolare, la garanzia non copre i danni causati da uso improprio, mancato rispetto delle istruzioni per l'uso, riparazioni e modifiche non autorizzate e usura.

APERA INSTRUMENTS (Europe) GmbH
Wilhelm-Muthmann-Straße 18, 42329 Wuppertal, Germany
Tel. +49 202 51988998 | info@aperainst.de | www.aperainst.de

